

Mask Aligner work protocol

1. All users must reserve the tool in BOOKIT.
2. Never turn on the front panel screen before turning on the controller.
3. Do not turn off the Mask aligner when your work is done (unless it is the end of the day and you sure that you are the last user on the Mask Aligner.
4. If the Mask Aligner is on "Flood exposure" mode and you are doing "load" your sample will be expose directly!
5. Always edit your parameters before loading your sample.
6. When your mask is on the chuck inside the Mask Aligner and you are inside "Change mask" mode never press on "Enter"-It will turn off the Mask vacuum and the mask will fall.