



1. PURPOSE

This Standard Operating Procedure (SOP) intends to describe the safe use of urethane in laboratory animals.

2. RESPONSIBILITY

Principal investigator (PI) and their research staff.

3. HAZARD CONSIDERATIONS

- 3.1. Carcinogen - Known to cause cancer in humans
 - 3.2. Irritant - eyes, mucous membranes, skin
 - 3.3. Toxicity - Kidneys, liver, CNS
-

4. PERSONAL PROTECTIVE EQUIPMENT (PPE)

- 4.1. Those handling urethane must wear chemically resistant gloves, a lab coat, eye protection, and appropriate lab attire (pants, closed-toe shoes).
 - 4.2. Urethane work should be done in a chemical fume hood.
 - 4.3. If manipulations occur outside of a chemical fume hood, personnel must wear a respirator equipped with a multi-gas filter until an exposure assessment can be conducted
-

5. PREPARATION OF URETHANE SOLUTIONS

- 5.1 Those preparing urethane solutions must always handle it inside a certified chemical fume hood.
 - 5.2 Care should not generate any aerosol during the preparation or injection procedure.
 - 5.3 Personal handling urethane must wear protective equipment (section 4.1)
-

6. ENVIRONMENTAL CONSIDERATION

- 6.1. Work with Urethane only in a chemical fume hood.
- 6.2. Care should be taken not to generate any aerosol during the preparation or injection.
- 6.3. Use a bench liner to contain the material.

- 6.4. Any leftover/unused urethane should be collected for disposal as chemical waste.
- 6.5. Used needles/syringes should be disposed of in a sharp's container destined for incineration, a cytotoxic sharps container.
- 6.5. Decontaminate the affected area with soap and water after absorbance in a spillover event.

7. ACCIDENT RESPONSE PROCEDURES

- 7.1. If the skin is exposed, wash immediately with soap and water. Flush mucus membranes with large amounts of water. Use a wet shower in case of extensive contamination.
 - 7.1.1. Report the incident to the supervisor.
 - 7.1.2. Reports the accident/injury to the Biosafety Unit.
 - 7.1.3. Remove all sources of ignition from the spill area.
 - 7.1.4. Spills in fume hood - use absorbent pads or vermiculite to clean up small fume hood spills— Clean up the spill area with additional pads or paper towels, followed by clean water.
 - 7.1.5. Spills in the room – notify others of the spill and keep the spill confined.
- 7.2. Spills must be cleaned immediately by adequately protected and trained staff.
- 7.3. Respiratory protection is required to clean up urethane spills outside an operating fume hood, as well as a disposable lab coat, goggles, and two pairs of nitrile gloves.
- 7.4. Use the same procedure as " spill in the fume hood."
- 7.5. Place spill waste in a plastic bag and label it as hazardous chemical waste.

8. REQUIRED AUTHORIZATION

- 8.1. The chemical safety officer must approve the suitability of the location and equipment to use urethane and have a startup meeting with the PI and research team before procurement.
- 8.2. The Principal Investigator (PI) must provide lab-specific training to all laboratory workers specific to the hazards (physical and health) involved in working with the substance, work area decontamination, and emergency procedures.
- 8.3. The PI must provide a copy of the MSDS and this SOP to any lab work with any materials covered.
- 8.4. The PI must ensure that all lab personnel has attended the required training and refresher training.

<p>SOP 601 URETHANE Revised on (dd,mm,yyyy) 01.11.2022 Approved by the BGU Animal Policy and Welfare Oversight Committee</p>
